pH 3310 IDS

DIGITAL METER FOR IDS SENSORS

a xylem brand

pH 3310 IDS - Contents

EN English	2
DE Deutsch.....	6
CS Český	6
DA Dansk	10
EL Ελληνικά	10
ES Español.....	12
FI Suomi	16
FR Français	16
HU Magyar	20
IT Italiano	20
JA 日本語	22
KO 한국어.....	26
NL Nederlands	26
NO Norsk	28
PL Polski.....	30
PT Português	32
RO Română	36
RU Русский	36
SK Slovenský	38
SV Svenska	40
TR Türkçe	42
ZH 中文.....	46
Contact Information	49

pH 3310 IDS

Overview

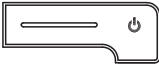
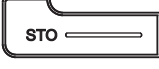
The pH 3310 IDS meter enables you to perform measurements (pH) quickly and reliably.

The pH 3310 IDS meter provides a maximum degree of operating comfort, reliability and measuring certainty.

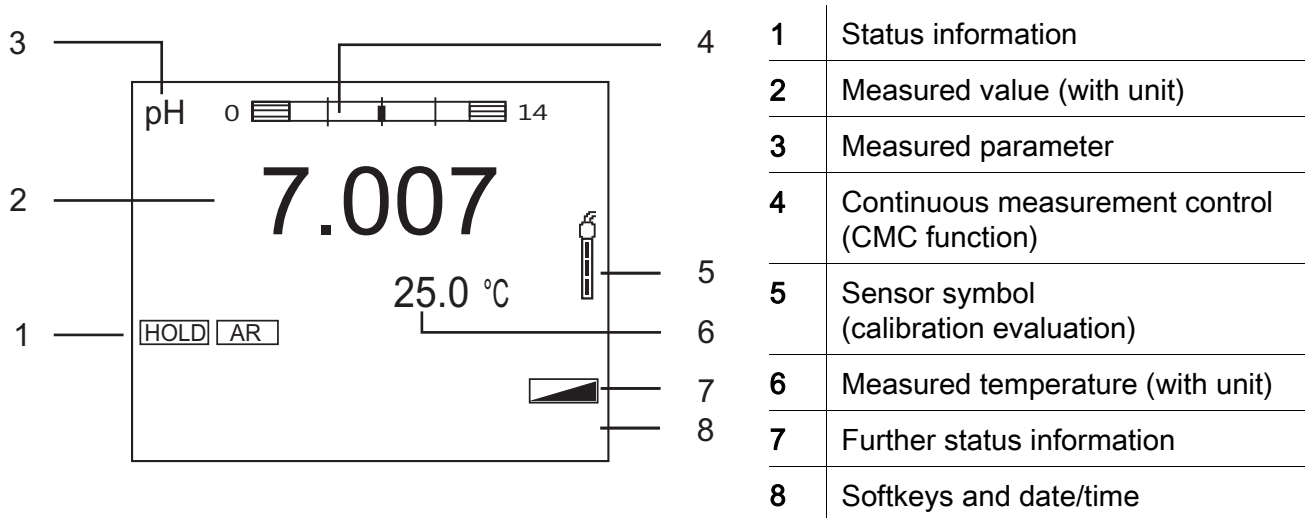
Keypad

Keys are indicated by brackets <.>.

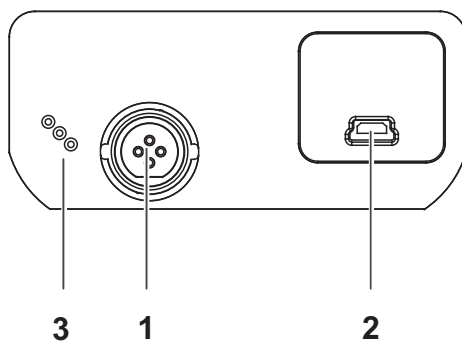
The key symbol (e.g. <ENTER>) generally indicates a short keystroke (under 2 sec). A long keystroke (approx. 2 sec) is indicated by the underscore behind the key symbol (e.g. <ENTER__>).

	<F1>:	Situation dependent function (is shown on the display)
	<F1__>:	<F1>/[Menu]: Opens the menu for measurement settings
	<F2>:	<F1__>/[Menu]: Opens the menu for system settings
	<F2__>:	
	<On/Off>:	Switches the meter on or off
	<M>:	Selects the measured parameter
	<CAL>:	Calls up the calibration procedure
	<CAL__>:	Displays the calibration data
	<STO>:	Saves a measured value manually
	<STO__>:	Opens the menu for the automatic save function
	<RCL>:	Displays the manually stored measured values
	<RCL__>:	Displays the automatically stored measured values
	<▲><▼>:	Menu control, navigation
		
	<ENTER>:	Opens the menu for measurement settings / confirms entries
	<ENTER__>:	Opens the menu for system settings
	<AR>:	Freezes the measured value (HOLD function) Switches the AutoRead measurement on or off

Display

Jack field /connections

1	pH electrode
2	Reference electrode
3	Temperature sensor
4	USB B interface (device)
5	Service interface

Initial commissioning

- 1 | Insert the supplied batteries. Make sure that the poles are in the correct position.
- 2 | Switch on the meter with the <On/Off> key.

Caution

Follow the safety instructions of the used sensors. The operating manuals of the sensors are available on the supplied CD and on the Internet under www.WTW.com.

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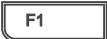
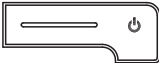
Display

The diagram shows the display interface with the following elements labeled:

- 1: Information about the current state (e.g., HOLD, AR)
- 2: Measured value (7.007) with unit
- 3: Parameter (pH)
- 4: Continuous control of measured values (CMC function)
- 5: Sensor symbol (evaluation of calibration)
- 6: Measured temperature (25.0 °C) with unit
- 7: More information about the current state
- 8: Softkeys and date / time

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The diagram shows the connection ports on the device:

- 1: pH Electrode
- 2: Reference Electrode
- 3: Temperature Sensor
- 4: USB-B Interface (device)
- 5: Service Interface

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The diagram shows a digital display with the following elements labeled 1 through 8:

- 1: HOLD and AR buttons
- 2: Large numerical pH reading (7.007)
- 3: pH scale bar (0 to 14)
- 4: CMC (Continuous Measurement Control) indicator
- 5: Sensor icon and calibration status
- 6: Temperature reading (25.0 °C)
- 7: Battery level indicator
- 8: Soft key and time/date display

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The diagram shows a socket panel with the following ports labeled 1 through 5:

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What can Xylem do for you?

We're a global team unified in a common purpose: creating innovative solutions to meet our world's water needs. Developing new technologies that will improve the way water is used, conserved, and re-used in the future is central to our work. We move, treat, analyze, and return water to the environment, and we help people use water efficiently, in their homes, buildings, factories and farms. In more than 150 countries, we have strong, long-standing relationships with customers who know us for our powerful combination of leading product brands and applications expertise, backed by a legacy of innovation.

For more information on how Xylem can help you, go to xylem.com.

®

zeller GmbH

Labworld.at Laborgeräte - Glas - Reagenzien
Mikrobiologie - Hygienekontrolle
Industriestr. 1, 6845 Hohenems, Austria
Tel. +43 (0)5576 76705 Fax +43 (0)5576 76705 7
Email: office@labworld.at